



Ractopamine ELISA Kit

*Competitive enzyme immunoassay kit for
quantitative analysis of Ractopamine*

Catalog Number: EL2052-01

For Research Use Only. Not for use in Diagnostic Procedures.

1. Background

Ractopamine (RAC) is a feed additive to promote muscle size in animals being raised as livestock for their meat (e.g. swine, cattle and turkeys) prior to their slaughter. It is in the beta-agonist family which has been associated with potential risk to human health. Therefore, Ractopamine has been banned for use as a livestock growth promoter in many countries around the world.

2. Test Principles

The Ractopamine plate kit is a competitive enzyme-labeled immunoassay. The Ractopamine sample extract and calibrators are pipetted into the test wells followed by the Ractopamine antibody into the test wells to initiate the reaction. During the 30 minutes incubation period, Ractopamine from the sample and Ractopamine antigen compete for binding to the Ractopamine antibody. The Ractopamine antibody is captured on the walls of the test well. Following this 30-minute incubation, the contents of the wells are removed and the wells are washed to remove any unbound Ractopamine and free Ractopamine antibody. After the wash step, IX HRP-conjugated Antibody#2 is added for 30 minutes incubation. The wells are washed afterwards, and a clear substrate is then added to the wells and any bound enzyme conjugate causes the conversion to a blue color. Following a 15-minute incubation, the reaction is stopped and the amount of color in each well is read. The color of the samples is compared to the color of the calibrators, and through this the Ractopamine concentration of the samples is derived.

3. Applications

This kit can be used for rapid testing of Ractopamine in urine, meat and feed.

4. Cross Reactions

Ractopamine	100%
Clenbuterol.....	< 0.1%
Salbutamol.....	< 0.1%

5. Equipment and Reagents Needed (not provided)

5.1 Equipment

- ELISA Reader (450nm/630nm)
- Deionized water
- Tissue Homogenizer
- Vortex mixer
- Timer
- Wash bottle
- Polystyrene centrifuge tube: 50mL, 2mL

- Micropipettes: 20 μ L-200 μ L, 100 μ L-1000 μ L
- 300 μ L-multipipette

6. Components Provided in This Kit

- Microtiter plate with 96 wells coated with Ractopamine
- Ractopamine Standards (6 vials \times 0.8mL/vial) 0ppb (green cap), 0.03ppb (purple), 0.1ppb (yellow cap), 0.3ppb (blue cap), 1ppb (orange cap), 5ppb (red cap)
- Ractopamine Antibody#1: 11mL
- 100X HRP-Conjugated Antibody#2: 0.25mL
- Antibody#2 Diluent: 20mL
- 20X Wash Solution: 28mL
- TMB Substrate Solution: 12mL
- Stop Solution: 14mL

7. Reagents Preparation

- 1X Wash solution: combine one volume of the 20X Wash Solution with 19 volumes of deionized water. Mix well.
- 1X HRP-conjugated Antibody#2: combine one volume of the 100X HRP-Conjugated Antibody#2 with 99 volumes of Antibody#2 Diluent. Vortex for 10 seconds to mix.
☛ Prepare this solution fresh before each test.

8. Notice and Precautions Before Operation

- Please use a fresh tip in the process of experiment and change the tips when absorbing different reagent.
- If running more than two strips at once, the use of a multichannel pipette is required.
- Make sure that all experimental instruments are clean.
- Treated samples can be stored at 2-8 $^{\circ}$ C for 24 hours in the dark.

9. Assay Process

9.1 Instructions Prior to Beginning Assay

1. Ensure that all reagents and microwells are at room temperature (20-25 $^{\circ}$ C).
2. Return all reagents to 2-8 $^{\circ}$ C immediately after their use.
3. Wash the microwells correctly; this is a vital factor in the reproducibility of the ELISA analysis.
4. Avoid direct sunlight during the incubation.

9.2 Steps in the Assay Process

1. Take all reagents out at room temperature (20-25°C) for more than 30 minutes. Shake gently before use.
2. Get the microwells needed out and return the rest into the zip-lock bag at 2-8°C immediately.
3. The diluted wash solution should be brought to room temperature before use.
4. Number every microwell position and all standards and samples should be run in duplicate. Record the standards and samples positions.
5. Dispense **50µL of Ractopamine Standards, positive control, or sample** into each well.
6. Dispense **100µL of Antibody#1** into appropriate test wells.
7. Shake the plate gently for 30 seconds using a back-and-forth motion.
8. Cover the plate. Incubate for **30 minutes** at room temperature.
9. Decant the contents of the wells into an appropriate waste container.
10. Rinse the microwells with 250µL of the 1X Wash Solution 3 times.
11. Absorb the residual solution by inverting with absorbent paper to remove the last of the Wash Solution.
12. Add **150µL of 1X HRP-Conjugated Antibody#2** (freshly prepared) to each well.
13. Shake the plate gently for 30 seconds using a back-and-forth motion.
14. Cover the plate. Incubate for **30 minutes** at room temperature.
15. Decant the contents of the wells into an appropriate waste container.
16. Rinse the microwells with 250µL of the 1X Wash Solution 3 times.
17. Absorb the residual solution by inverting with absorbent paper to remove the last of the Wash Solution.
18. Add **100µL TMB Substrate Solution** to each well, mix gently by shaking the plate manually and incubate for **15 minutes** at 25°C with cover.
19. Add **100µL of Stop Solution** to each well. Mix gently by shaking the plate manually and measure the absorbance at 450nm (Read the result within 5 minutes after addition of Stop Solution).

10. Results

10.1 Calculating the Percentage absorbance

- The mean values of the absorbance values obtained from the standards and the samples are divided by the absorbance value of the first standard (zero standard) and multiplied by 100%.

$$\text{Absorbance (\%)} = B / B_0 * 100$$

B = the mean absorbance value of each standard or each sample

B₀ = absorbance value of zero standard

10.2 Drawing a Standard Curve

- To draw a standard curve, the absorbance value of standards as y-axis, semilogarithmic of the concentration of the standards (ppb) as x-axis.
- The concentration of each sample (ppb), which can be read from the standard curve, is multiplied by the corresponding dilution factor of each sample followed, and the actual concentration of sample is obtained.
- Sample dilution factor: If the absorbance of a sample is lower than the highest calibrator (5 ppb), the concentration of Ractopamine is too high and out of range of the standard

curve. Dilute the sample and rerun. Samples should be diluted to fit into the standard curve (0.03 ppb to 5 ppb). Results must then be multiplied by the dilution factor used.

11. Sensitivity, Accuracy and Precision

11.1 Test Sensitivity:

- Overall Sensitivity..... 0.03ppb

11.2 Detection limit:

- Feed 1ppb
- Meat 0.1ppb
- Urine 0.4ppb

11.3 Accuracy:

- Feed, meat and urine $80 \pm 10\%$

11.4 Precision:

- C.V. of the ELISA kit less than 10%

12. General Instructions

12.1 Temperature of Reagents and Samples

- The mean values of the absorbance values obtained for the standards and the samples will be reduced if the reagents and samples have not been restored to room temperature (20-25°C).

12.2 Microwells

- Do not allow Microwells to dry between steps to avoid unsuccessful reproducibility and operate the next step immediately after tapping the microwells holder.

12.3. Shaking of Reagents

- Shake each reagent gently before use.

12.4. Skin Protection

- The Stop Solution is 0.75N HCl, keep your skin away from it.

12.5 Out of Date Kits

- Don't use kits that are expired. Don't exchange the reagents of different batches, or else it will drop the sensitivity.

12.6 General Comments

- Keep the ELISA kits at 2-8°C, do not freeze. Store the unused microwell plates back to the foil pouch. Avoid straight sunlight during all incubations. Covering the microtiter plates is recommended.

12.7 Special Issues Concerning Solutions and Reagents

- TMB Substrate should be discarded if it turns a blue hue, as it should be clear. One or more of the reagents may be bad if the absorbance value of the zero standard is less

than 0.5 ($A_{450\text{nm}} < 0.5$).

12.8 Special Issues Concerning Color

- The coloration reaction takes 15 minutes after the addition of TMB Substrate, but you can prolong the incubation time ranges to 35 minutes or more if the color is too light to be determined, never exceed 40 minutes, on the contrary, shorten the incubation time properly.

12.9 Incubation Temperatures

- Incubation temperature should be at room temperature (20-25°C). Higher or lower temperature on the day of testing will lead to experiment-to-experiment changes.

13. Storage

- Storage condition: 2-8°C
- Storage period: 12 months or until expiration date
 - For most accurate results, only use reagents and standards from the same kit. The kit's expiration date is thereby that of the earliest expiring component.

Customer Notes:

Contact Us

Attogene

3913 Todd Lane, Suite 310 Austin, TX 78744

Phone: 512-333-1330

Email: sales@attogene.com

Web: www.attogene.com

EL2052-01__v.2